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PSC_SC_700310T_IQ-HydroFlex

Purpose:

This form serves the TECAN authorized Service Technician as a checklist on the main installation process. However, this form does not contain the procedure in detail (refer to reference documents and to the SOP Field Intervention Doc. 10402TMs01). In the following sections place a checkmark in the check box that applies. Actions flagged N/A must be explained in the notes section of this document if the reason differs from the description text. Skipped or failed actions must always be explained. Printed versions of this checklist must always be bundled with the corresponding associated documents.

Disclaimer:

Depending on the configuration of instrument, sections of checklist can be omitted and may not be displayed.

1. General Information

Document Title IQ-HydroFlex
 Document Number SC 700310T
 Document Version 06

Customer Information

Service Order # 000004507113
 Contact person PROSPERI ALICE
 Customer e-mail alice.prosperi@izsler.it
 Performed By Salvatore Trolano (salvatore.trolano@tecan.com)
 Lab site Laboratorio P2 01 Floor 2

Instrument Information

Instrument I.Z.S.Lombardia e Emilia Romagna; INSTRUMENT HYDROFLEX PLUS
 Code 108401-161937; 12481170
 Serial Number ---; 2303011733
 Firmware Version V 4.22
 Instrument Installation New Installation

2. Configuration

Vacuum Filtration NO
 Channel 4
 MBS NO
 LLD NO
 HydroControl NO

3. Reference documents

Document Instructions for Use HydroFlex

4. Site check and delivery control

Site check according to specifications, see instructions for use ☒ YES ☐ NO
 (IFU) HydroFlex
 All items delivered and functional according to IFU HydroFlex ☒ YES ☐ NO
 and IFU HydroFlex Options

5. Installation of Instrument

Visual check for transport damage ☒ YES ☐ NO
 Unpacking & positioning of instrument ☒ YES ☐ NO
 Installation of the instrument according to IFU HydroFlex ☒ YES ☐ NO

6. Installation of Liquid Connections

Connect the waste system, see IFU HydroFlex ☒ YES ☐ NO
 Connect the solution tubes, see IFU HydroFlex ☒ YES ☐ NO

8. Customer instruction

Instruction of Instrument Warning Labels Done
 How to connect Washer with Software Done
 Basic Explanation Software Menu Structure according to IFU Done
 Maintenance -- according to the manufacturer's instruction to be performed every 12 months Done
 List of Instructed Employees at customer company

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11. Signatures

Confirmation

The service technician confirms with signature that the intervention was performed in accordance to this checklist and the published Tecan procedures that apply to the instrument listed on this form.

Date

02.05.2023

Checklist specific signature is NOT required

☒ YES

☐ NO

Service Bundle

☒ YES

☐ NO

Doc. No.	Title	Version	Effective Date	Author
SC_700310T	IQ-HydroFlex	06	2021-07-19	Markus Relschenböck

© Copyright Tecan Austria GmbH • Unterebergstr. 1a • 5082 GRÖDIG / AUSTRIA

PSC_SC_700330T_OQ-HydroFlex

Purpose:

This form serves the TECAN authorized Service Technician as a checklist on the main verification process. However, this form does not contain the procedure in detail (refer to reference documents and to the SOP Field Intervention Doc 10402TMs01). This checklist serves for hardware verification only and does not cover any process validation. Actions flagged N/A must be explained in the notes section of this document if the reason differs from the description text. Skipped or failed actions must always be explained. Printed versions of this checklist must always be bundled with the corresponding associated documents.

Disclaimer:

Depending on the configuration of Instrument, sections of checklist can be omitted and may not be displayed.

1. General Information

Document Title	OQ-HydroFlex
Document Number	SC_700330T
Document Version	08

Customer Information

Service Order #	000004507113
Contact person	PROSPER ALICE
Customer e-mail	alice.prosper@izsler.it
Performed By	Salvatore Trolano (salvatore.trolano@tecan.com)
Training Certificate attached	☒ YES ☐ NO
Lab site	Laboratorio P2 01 Floor 2

Instrument Information

Instrument	I.Z.S.Lombardia e Emilia Romagna; INSTRUMENT HYDROFLEX PLUS
Code	108401-161937; 12481170
Serial Number	—; 2303011733
Balance Serial No.	B909318773
Calibration Certificate attached	☒ YES ☐ NO
Service Software Version	V 4.5.4
Firmware Version	V 4.22
Instrument Installation	New installation

2. Configuration

Vacuum Filtration	NO
Channel	4
MBS	NO
HydroControl	NO

3. Reference documents

Document	Instructions for Use HydroFlex
----------	--------------------------------

4. Operation Values

Proceed to the next chapter.

4.1. Installation Qualification (IQ)
Check if IQ is available

Acceptance values	IQ is available and all applicable test points are "done" or "passed"
Status	Passed

4.2. Function Test Channel 1

Define the following program to perform a function test:

1. Plate Mode
2. Plate Type: Greiner 96-well flat bottom
3. One Cycle
4. One Asp. step with the following parameters: Mode: NORMAL; Pos: BOTTOM; Time: 1 s; Head speed: 10 mm/s; Asp. Rate: 2
5. One Disp. step with the following parameters: Pos: OVERFLOW; Volume: 300 µl; Channel: 1; Disp. Rate: 300 µl/s
6. One Soak step with the following parameters: Time: 1:30 min; Shake: YES; Intensity: MEDIUM
7. One Wash step with the following parameters: Pos: OVERFLOW; Vol: 3000 µl; Channel: 1; Head speed: 10 mm/s; Wash Rate: 300 µl/s
8. Final aspirate with the following parameters: Mode: CROSSWISE; Pos: BOTTOM; Time: 4 s; Head speed: 10 mm/s

Acceptance values	Function test if finished without errors.
Status	Passed

4.3. Function Test Channel 2

Define the following program to perform a function test:

1. Plate Mode

2. Plate Type: Greiner 96-well flat bottom

3. One Cycle

4. One Asp. step with the following parameters: Mode: NORMAL; Pos: BOTTOM; Time: 1 s; Head speed: 10 mm/s; Asp. Rate: 2

5. One Disp. step with the following parameters: Pos: OVERFLOW; Volume: 300 µl; Channel: 2; Disp. Rate: 300 µl/s

6. Final aspirate with the following parameters: Mode: CROSSWISE; Pos: BOTTOM; Time: 4 s; Head speed: 10 mm/s

Acceptance values

Function test if finished without errors.

Status

Passed

4.4. Function Test Channel 3

Repeat the "Function Test Channel 2" using Channel 3 in the dispense and wash steps

Acceptance values

Function test is finished without errors.

Status

Passed

4.5. Function Test Channel 4

Repeat the "Function Test Channel 2" using Channel 4 in the dispense and wash steps

Acceptance values

Function test is finished without errors.

Status

Passed

4.8. Dispensing Accuracy

Follow the instructions for "Dispense Accuracy Check" in the Instructions for Use HydroFlex (Chapter Performance Testing / Quality Control)

Acceptance values

Weight of dispensed liquid must be ≥ 18.8 g and ≤ 19.8 g.

Measured Value

19.52

Status

Passed

4.9. Residual Volume

Follow the instructions for "Residual Volume Check" in the Instructions for Use HydroFlex (Chapter Performance Testing / Quality Control)

Acceptance values

Weight of dispensed liquid must be ≤ 0.190 g.

Measured Value

0.13

Status

Passed

6. Signatures

Confirmation

The service technician confirms with signature that the intervention was performed in accordance to this checklist and the published Tecan procedures that apply to the instrument listed on this form.

Date

02.05.2023

Due date next maintenance

02.05.2024

Checklist specific signature is NOT required

☒ YES

☐ NO

Service Bundle

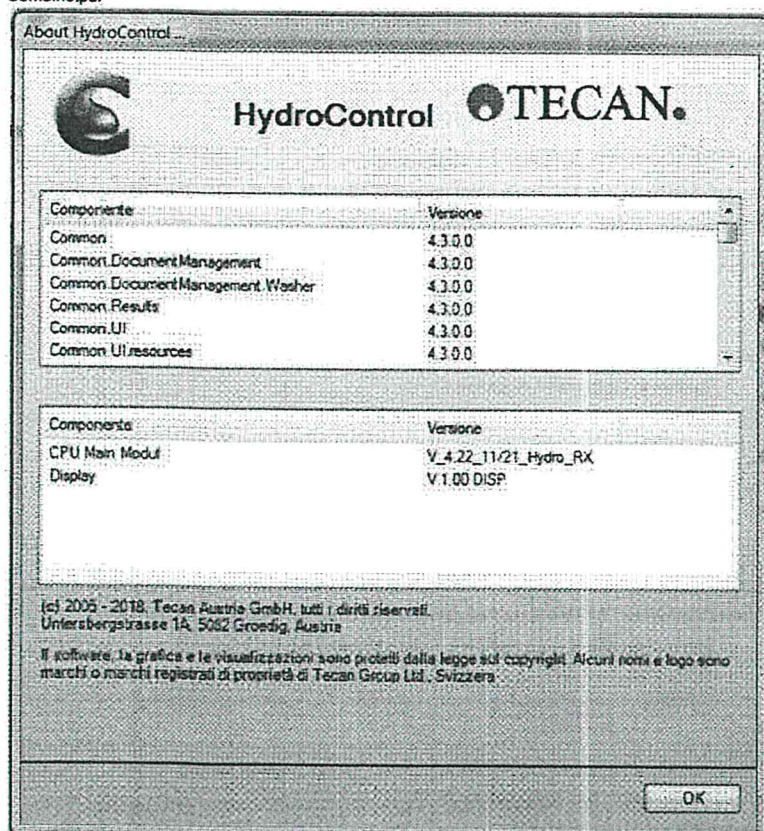
☒ YES

☐ NO

Doc. No.	Title	Version	Effective Date	Author
SC_700330T	OQ-HydroFlex	08	2021-07-19	Markus Reischenböck

© Copyright Tecan Austria GmbH

Combine.pdf

Certificato di Taratura N°
IT0069-040-072122-ACC

METTLER TOLEDO

Mettler Toledo S.p.A.
Via Anna Maria Mozzoni, 2/1
I-20152 Milano, Italy
Nr. Verde 800300693
assistenza@mt.com

Accuracy Calibration Certificate

Accuracy Calibration Certificate

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Tecan Service Report 4507113 for I.Z.S. Lombarde e Emilia Romagna.pdf

Certificato di Taratura N°
IT0089-040-072122-ACC

METTLER TOLEDO Service

Risultati

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
Test Load: 10 g

	As Found	As Left
1	9,99987 g	10,00002 g
2	9,99986 g	10,00003 g
3	9,99988 g	10,00002 g
4	9,99988 g	10,00004 g
5	9,99989 g	10,00003 g
6	9,99985 g	10,00004 g
7	9,99986 g	10,00004 g
8	9,99986 g	10,00004 g
9	9,99987 g	10,00004 g
10	9,99987 g	10,00004 g

Deviazione Standard	0,000012 g	0,000008 g
---------------------	------------	------------

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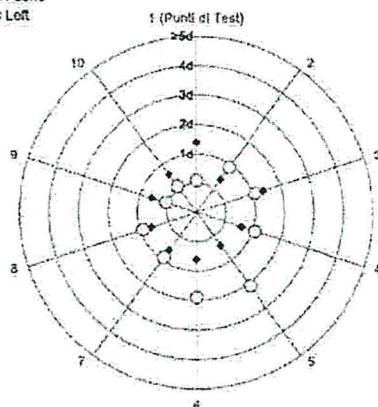
Test Load: 100 g

Posizione	As Found	As Left
1	99,99968 g	99,99968 g
2	99,99963 g	99,99968 g
3	99,99962 g	99,99967 g
4	99,99959 g	99,99968 g
5	99,99961 g	99,99969 g

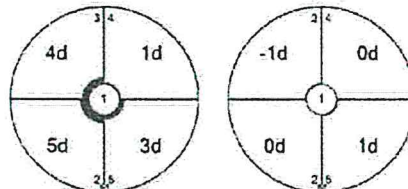Dev. Massima	0,00005 g	0,00001 g
--------------	-----------	-----------

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